


ARCspectro FT-MIR

AR@ptix
Switzerland

DATA SHEET



Compact fibered MIR 2.0-6.0 Fourier-transform spectrometer

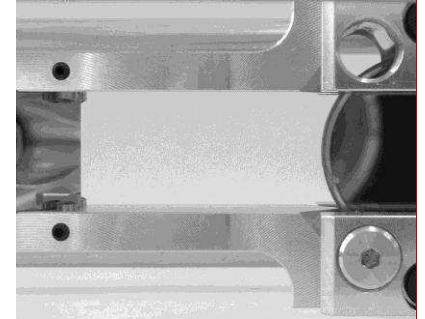

The FT-MIR "Rocket" is a fibered MIR spectrometer (2-6 microns wavelength) that is easily coupled to accessories using mid-infrared CIR fibers. It incorporates Arcoptix's 3rd generation new self-compensated interferometer, using a solid-state internal reference laser. The system features excellent sensitivity using a 2-TE cooled MCT detector, stability and wavelength reproducibility.

Typical characteristics of the FT-MIR series are :

- **Compact (world smallest FT-MIR)**
- **Broad spectral range 5000-1600cm⁻¹ (2.0–6.0 μm)**
- **Fiber input (SMA-905)**
- **High resolution of 4cm⁻¹**
- **Cost-effective**

Typical applications of the FT-MIR series are :

- **Chemometrics**
- **Material identification**
- **Laser characterization**
- **Gas analysis**

For additional information please contact:

info@taisl.com

<http://www.taisl.com/>

FEATURES & BENEFITS


Extreme wavelength range, high resolution and optical throughput are combined in our compact Fourier-Transform spectrometers

High performance interferometer

The ARCSpectro FT-MIR uses a compact yet high-performance, optically self-compensated interferometer.

Solid-state control laser

The use of a solid state reference laser allows our FT-MIR to be cost-effective, while maintaining excellent stability and wavelength reproducibility.

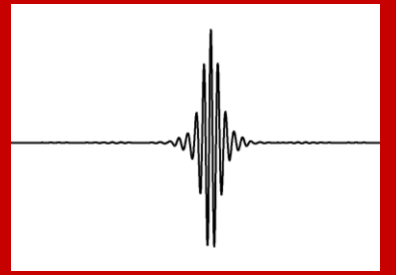


ARCSpectro FT-MIR

AR@ptix
Switzerland

DATA SHEET

Type of interferometer	Self-compensated multi-pass reciprocating interferometer
Spectral range	5000–1600cm ⁻¹ (2.0–6.0 μm)
Detector	2-stage TE-cooled MCT (HgCdTe) detector
Resolution	4 cm ⁻¹
Dynamic range	>1000:1
Measurement time	4 s (including processing)
Control laser	Solid-state 850nm
Wave-number accuracy	<1cm ⁻¹ @ 4000cm ⁻¹ (25± 5 °C)
Optical fiber input	SMA 905, fiber core diameter 500um, NA=0.2
Operating temperature	10 - 40°C
ADC resolution	24 bits
Power requirements	15 V / 1A (line-powered)
Communication interface	USB 2.0
Included software interfaces	Windows XP/Vista/7 software
Size	180mm x 126mm x 78mm
Weight	1700 g

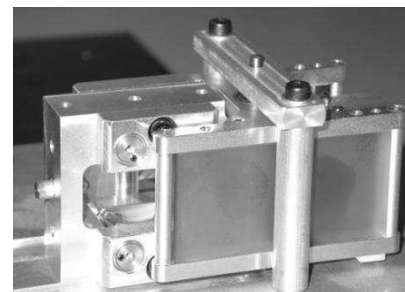

OPTIONS

Temperature-stabilized control laser
for ultimate wave-number accuracy down to <0.1cm⁻¹

External detector module
Alternative configuration, where the MIR light source is integrated into the interferometer, and an external detector is used.

Detector s
Other detectors are available on request.

Advanced software
Software modules are available for:

- Library search
- Chemometrics

Técnicas Asociadas de Instrumentación, S.L..

For more information about ARCOptix, visit <http://www.taisl.com/> Tel / Fax: +34 (91) 3555802

Specifications are subject to change without notice

© 2011 ARCOptix S.A.